




Products for Anatomic Pathology

### Procedure Super Pap Pen Liquid Blocker

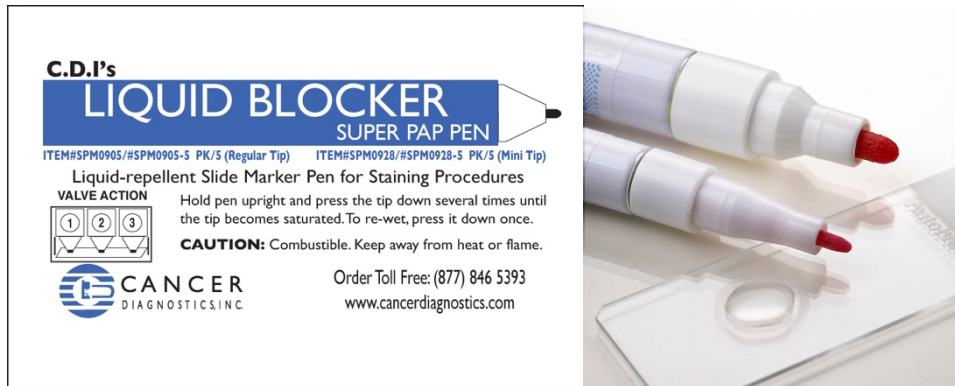

Before using pen/applying to slide, hold the pen point up and press it down with fingertip to remove excess air out of pen barrel to prevent overflow.

We recommend use on non-coated/charged slides.

If the glass slide is hydrophile treated, (coated/charged) the glass slide should be soaked in a 0.1N hydrochloric acid solution for about 30 seconds and then dried after washing with water to avoid “fall-off” of the circle from the slide.

#### For Frozen Sections:

Before applying the pen circle, it is necessary to clean the glass slide by xylene and ethanol or coat with poly-L-lysine.